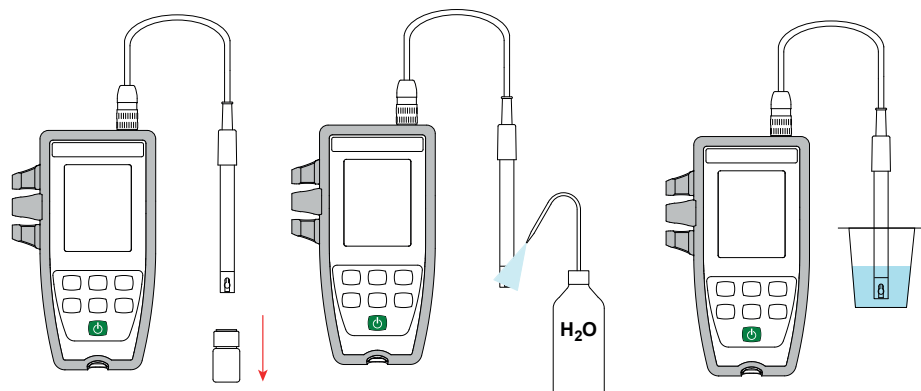
MESURE

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

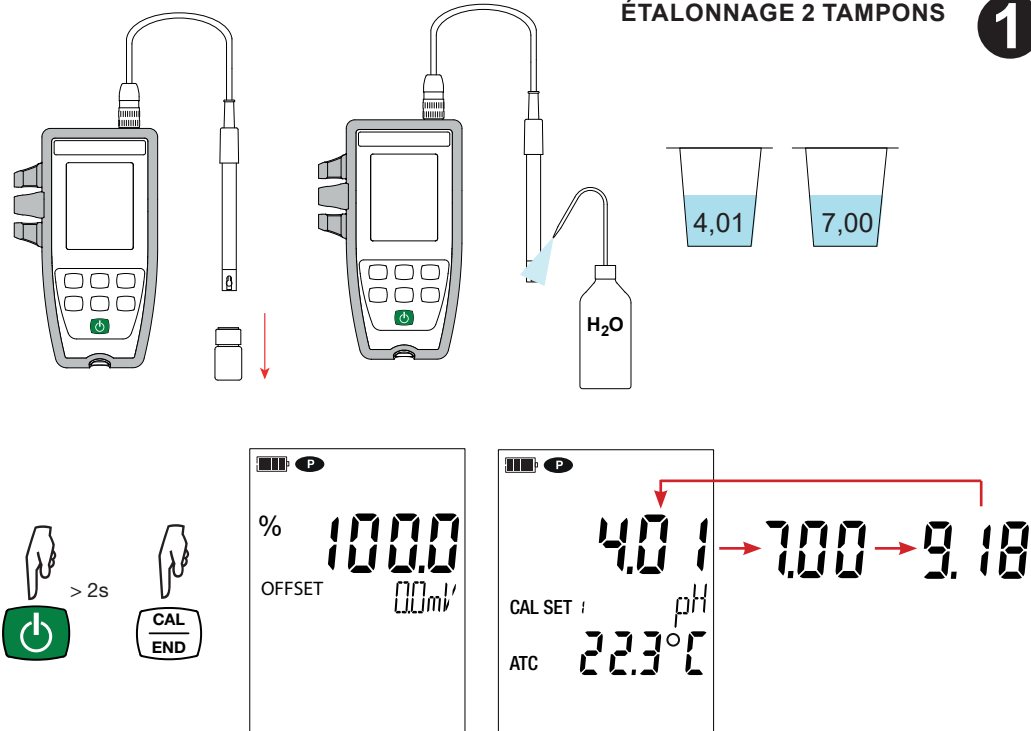
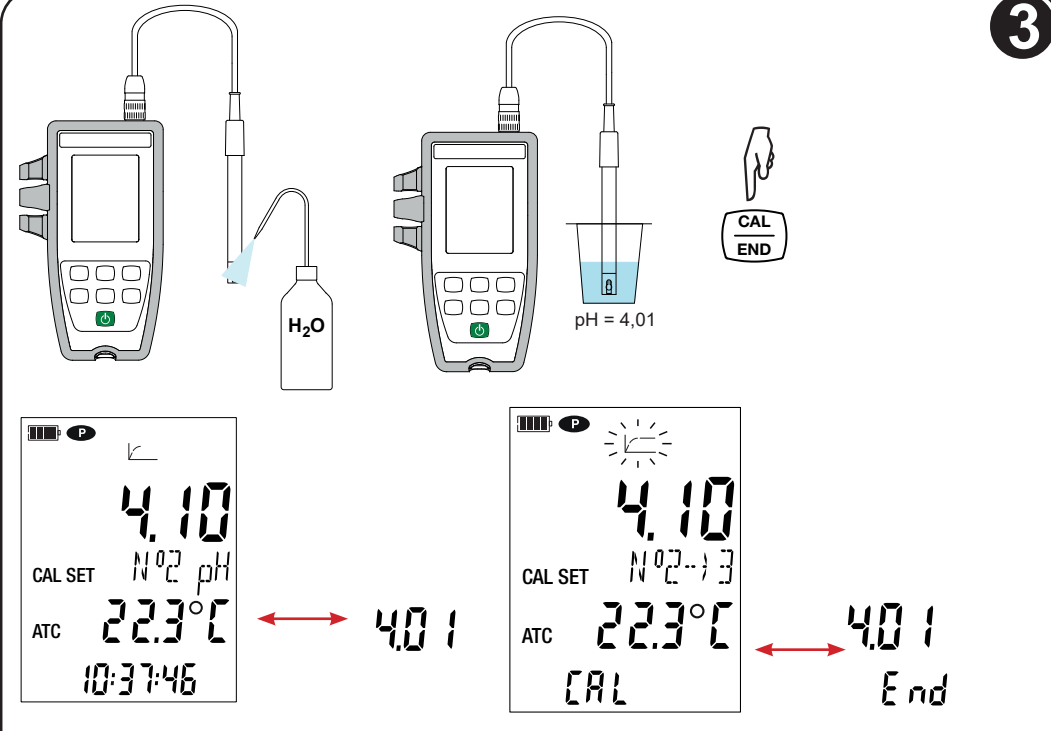
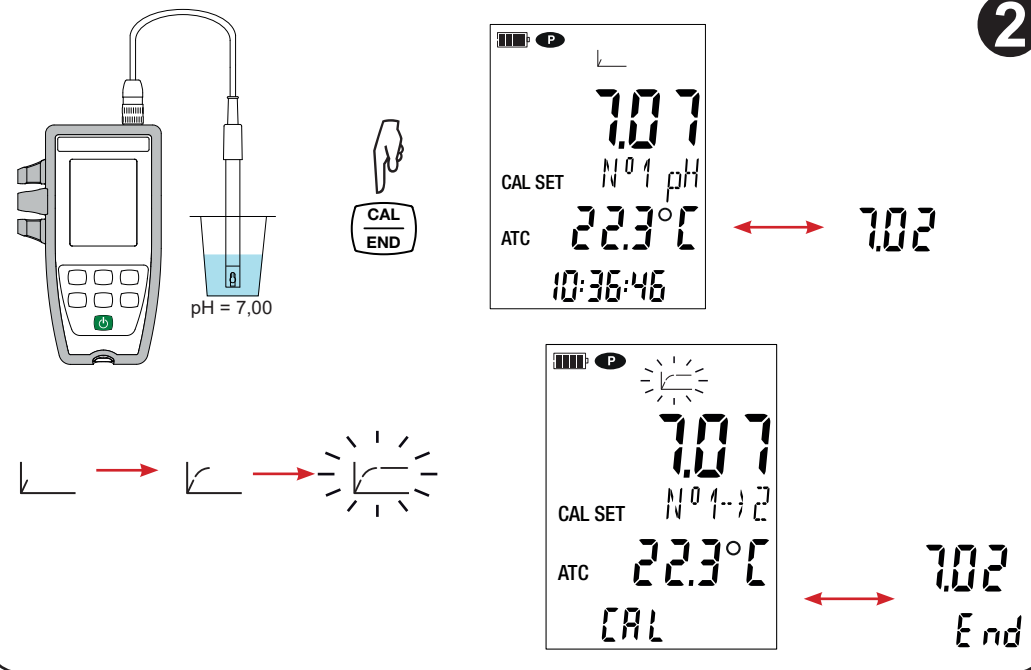
4

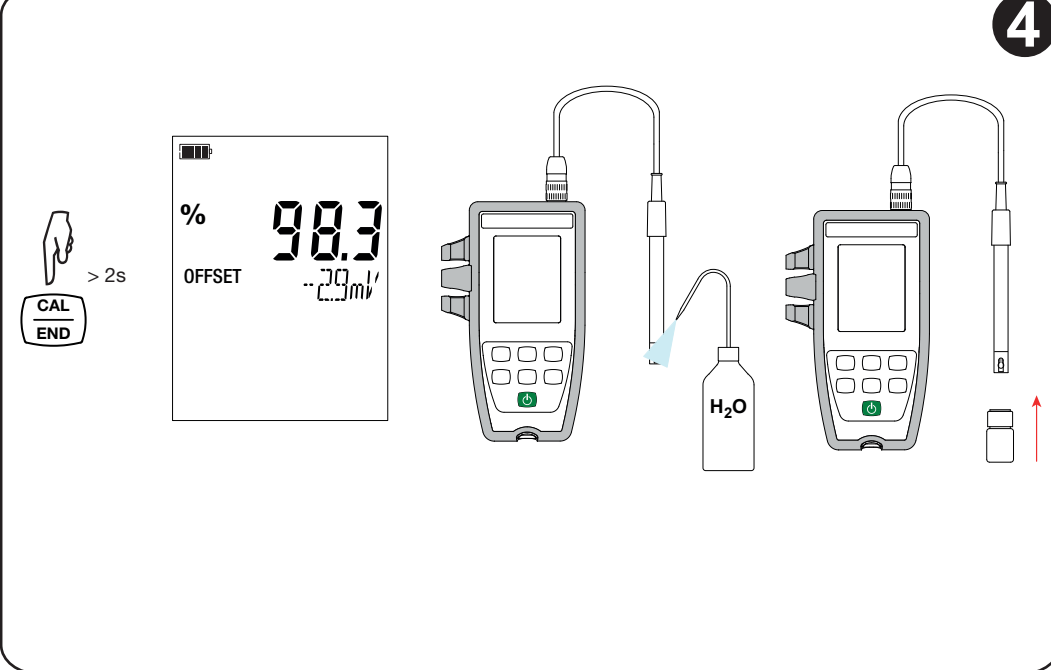MEASUREMENT

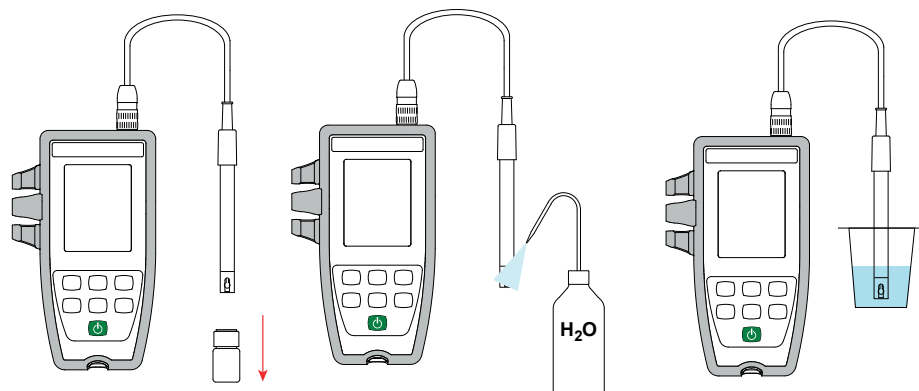The  symbol indicates the quality of the last electrode calibrated.

Slope value	Offset value	State of the electrode
$95\% \leq \text{slope} \leq 105\%$	offset ≤ 19 mV	
$90\% \leq \text{slope} < 95\%$	$19 \text{ mV} < \text{offset} \leq 38 \text{ mV}$	
$85\% \leq \text{slope} < 90\%$	$38 \text{ mV} < \text{offset} \leq 58 \text{ mV}$	
slope $< 85\%$ or slope $> 105\%$	$58 \text{ mV} < \text{offset}$	

FRANCE
Chauvin Arnoux
12-16 rue Sarah Bernhardt
92600 Asnières-sur-Seine
Tél : +33 1 44 85 44 85
info@chauvin-arnoux.com
www.chauvin-arnoux.com

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Measure up

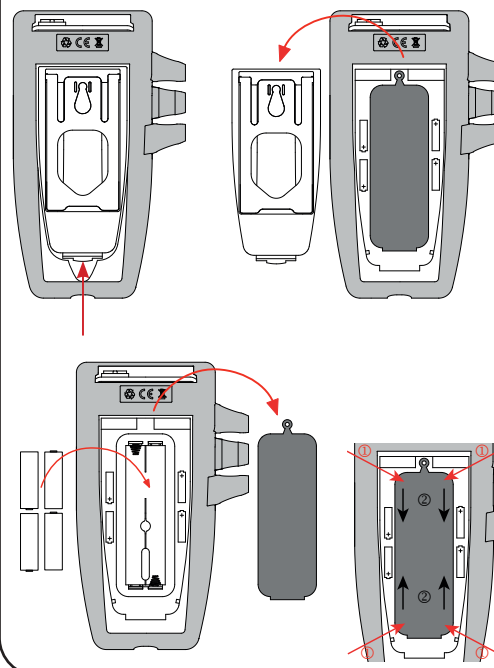
USER'S MANUAL

Visit our web site to download the operating instructions for your instrument and your electrodes:
www.chauvin-arnoux.com

Go to the **Products** tab and search on the name of your instrument. When you have found the instrument, go to its page. The operating instructions are on the right. Download them.

Proceed in the same way for the pH and ORP electrodes.

INSERTING THE BATTERIES

DATA LOGGER TRANSFER SOFTWARE

Visit our web site to download the latest version of the application software:
www.chauvin-arnoux.com

Go to the **Support** tab, then **Download our software**. Then search on the name of your instrument.

Download the software, then install it on your PC.

KEYS

Key	Function
	On/Off key (long press).
MEM REC	MEM: a short press to record the measurement and the temperature. REC: a long press to start or stop recording.
pH/mV °C/°F	pH/mV: a short press to select the type of measurement (pH or ORP). °C/°F: a long press to select the unit of temperature.
	Backlight.
CAL END	CAL: a short press to look up the calibration parameters and start a calibration. END: a long press to terminate the calibration.
 	- A short press to set the temperature in MTC or the parameters in SET mode. - A long press to select the standard solution during the calibration.

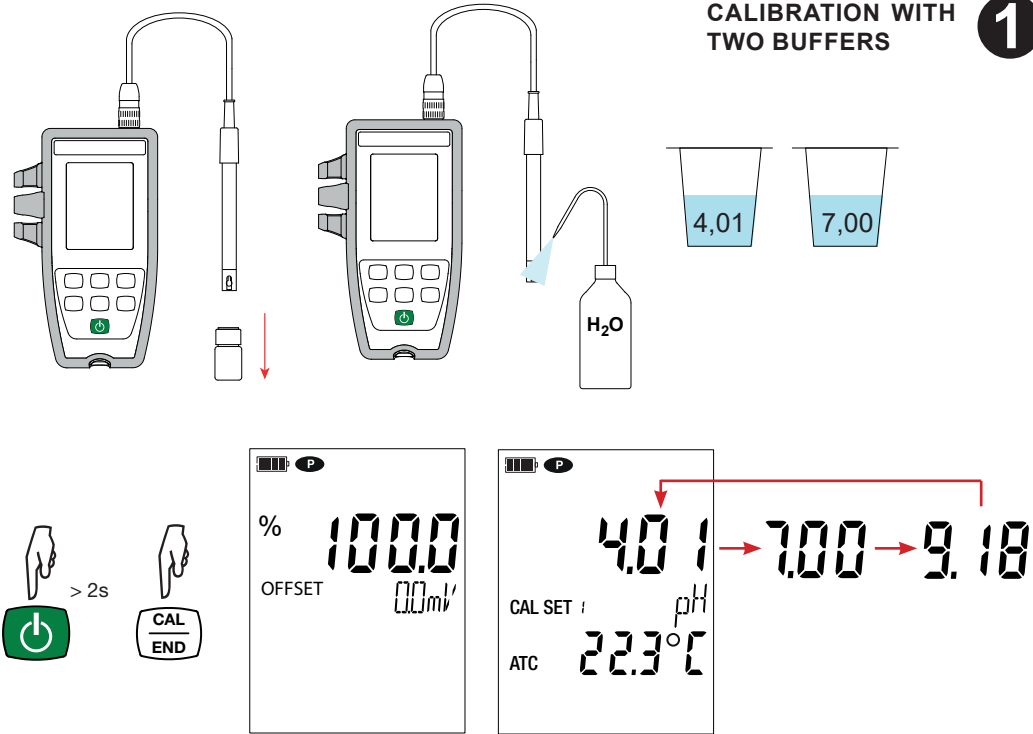
DELIVERY CONDITION

The C.A 10101 pH-meter is delivered in a carrying case with:

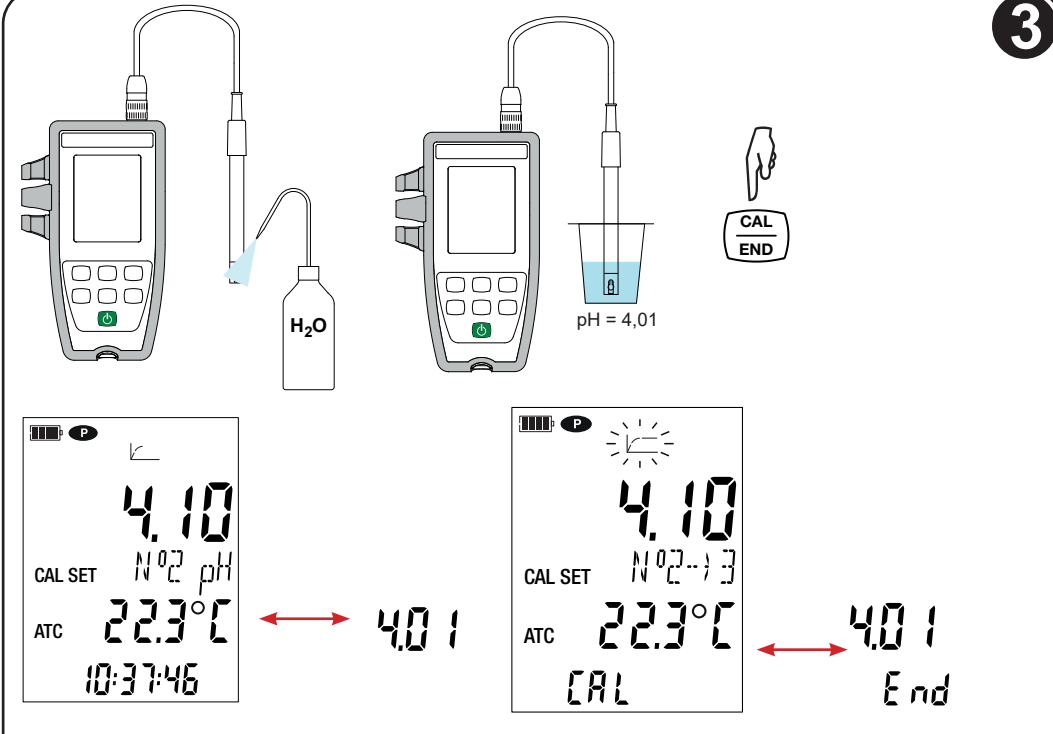
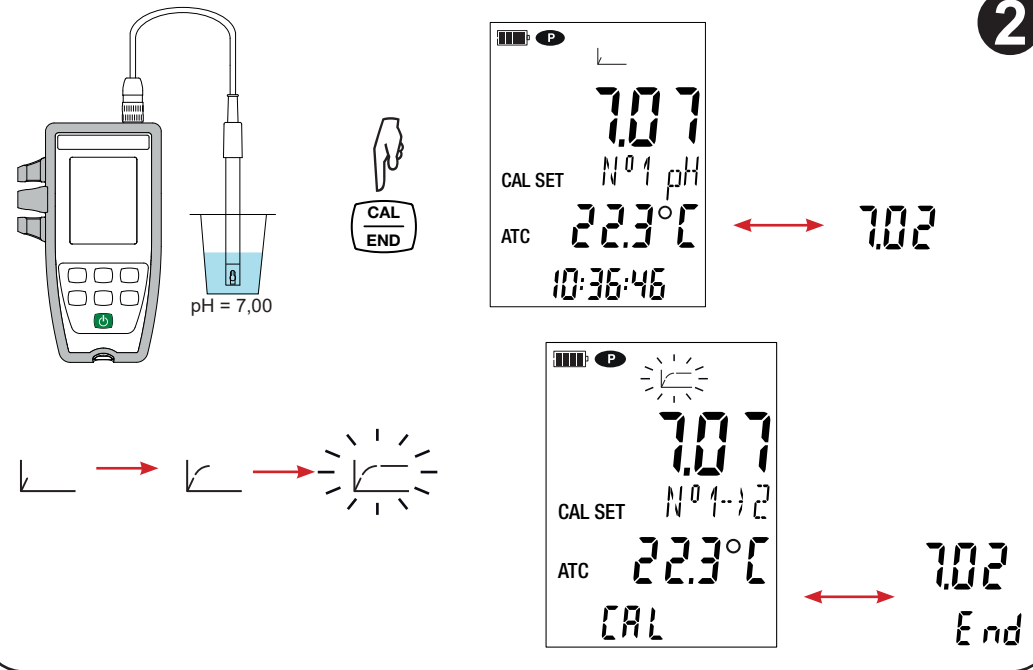
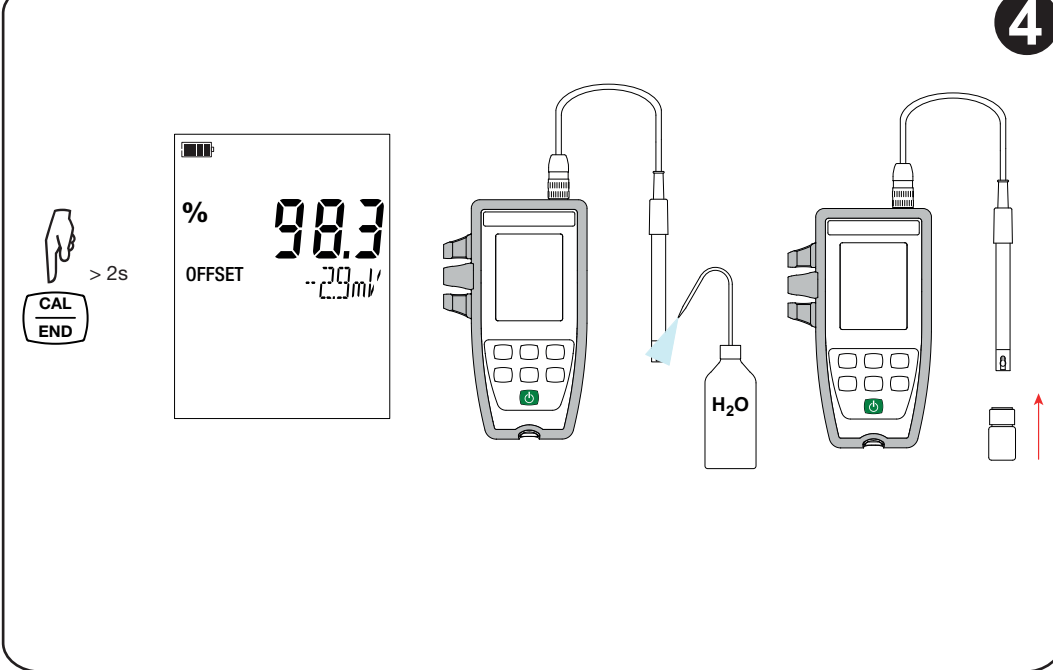
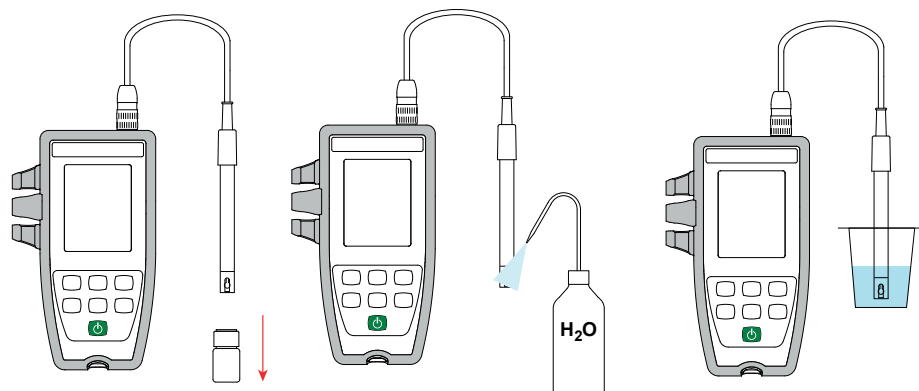
- One XRGST1 combined pH electrode with integrated temperature probe and storage reservoir.
- Four AA or LR6 alkaline batteries.
- One protective sheath fitted on the instrument.
- Two ready-to-use NIST buffer solutions, pH 4.01 and 7.00.
- Two beakers.
- One USB to micro USB cable.
- One wrist strap.
- Getting started guides (one per language).
- One test report.

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